




QC for Cell-Loaded HIVES

Receiving Instructions

- On dry ice, cell-loaded HIVES should be checked to make sure:
 - Dry ice remains in the package, any temperature sensors have not been activated, and the HIVE is still frozen
 - The liquid in the HIVE is completely covering the HIVE array and is white in color
 - The clear stopper is still present
- Cell-loaded HIVES should be placed at -80 immediately to ensure they remain frozen

Before Processing

- After HIVES are thawed for processing, check to make sure:
 - HIVES are not leaking
 - Liquid completely covers the array with no dry spots
 - Liquid is not present in the outside ring of the HIVE collector

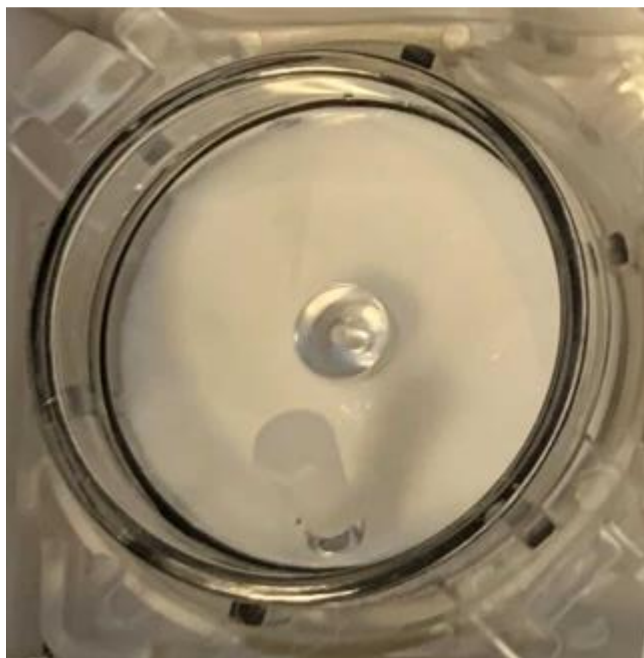

Good

(Liquid is frozen and completely covers the array)